

BrightStar™ BioDetect™ (Cat. # 1930)
Instruction Manual

- I. Introduction 1**
 - A. Overview
 - B. Reagents Provided with the Kit and Storage
 - C. Materials Not Provided with the Kit
 - D. Related Products Available from Ambion
- II. BioDetect Protocol 3**
 - A. When to Use BioDetect
 - B. Buffer and Equipment Preparation
 - C. Detection Procedure
- III. Troubleshooting 6**
 - A. Use of the Labeled Control DNA
 - B. High Background
 - C. Low Sensitivity
- IV. Additional Procedures 10**
 - A. Electroblothing from Polyacrylamide Gels
 - B. Crosslinking
- V. Appendix 13**
 - A. Specification Sheet
 - B. IX Blocking Buffer Material Safety Data Sheet
 - C. SDS Material Safety Data Sheet

Manual Version 0105

Literature Citation

We would appreciate if when you are describing a procedure using this product in a Materials and Methods Section for publication that you refer to it as the BrightStar™ BioDetect™ Kit.

If a paper that cites one of Ambion's products is published in a research journal, the author(s) may receive a free Ambion T-shirt by sending in the completed form at the back of this instruction manual, along with a copy of the paper.

Warranty and Liability

Ambion is committed to providing the highest quality reagents at competitive prices. Ambion warrants that the products meet or exceed the performance standards described in the product specification sheets. If you are not completely satisfied with any product, our policy is to replace the product or credit the full purchase price and delivery charge. No other warranties of any kind, expressed or implied are provided by Ambion. Ambion's liability shall not exceed the purchase price of the product. Ambion shall have no liability for direct, indirect, consequential or incidental damages arising from the use, results of use, or inability to use its products. This product is intended for research use only. This product is not intended for diagnostic or drug purposes.

Copyright pending by Ambion, Inc. all rights reserved.

CDP-Star® is a registered trademark of Tropix and is manufactured under U.S. Patent #5,330,900.

I. Introduction

A. Overview

The BrightStar™ BioDetect™ Kit for nonisotopic detection of biotinylated RNA and DNA probes is a result of research efforts at Ambion to develop a nonisotopic chemiluminescent detection system that does not suffer from the high background and low sensitivity typical of these systems. The result is a kit that yields extremely high signal to noise ratios with a limit of detection of about 100 fg.

The kit is designed to detect biotinylated RNA or DNA probes on positively charged nylon membrane. Ambion's BrightStar-Plus™ positively charged nylon membrane will give optimal performance with this chemiluminescent detection system.

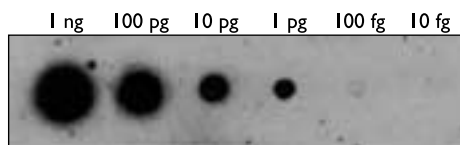


Figure 1. Detection of biotinylated DNA with BioDetect

A dilution series of a 1 kb biotinylated DNA probe, as indicated in the figure, was spotted on Ambion's BrightStar-Plus™ nylon membrane and then detected using the BrightStar™ BioDetect™ protocol. A 45 minute exposure to film revealed a limit of sensitivity of about 100 fg.

B. Reagents Provided with the Kit and Storage

Amount	Component	Storage
1 L	5X Wash Buffer	RT
3 L	1X Blocking Buffer	RT
125 ml	10X Assay Buffer	4°C
20 µl	Streptavidin-Alkaline Phosphatase*	4°C
50 ml	CDP-Star®†	RT
20 µl	Labeled Control DNA (1 ng/µl)*	-20°C

* These components are shipped in the slit of the grey foam inside the styro-foam cooler.

† CDP-Star® is a registered trademark of Tropix and is manufactured under U.S. Patent #5,330,900.

**IMPORTANT!**

Do not freeze the Streptavidin-Alkaline Phosphatase. Repeated freezing and thawing can compromise its performance.

Properly stored components are guaranteed for 6 months from the date received.

C. Materials Not Provided with the Kit

- Positively charged nylon membrane
We recommend Ambion's BrightStar™-Plus Membrane (Cat #10100, 10102, 10104)
- Electrotransfer device
Although virtually any electroblotter should function well, we recommend the Panther™ Semi-Dry Electroblotter (Model #HEP-1) from OWL Scientific as an affordable device that has been tested at Ambion for this application.
- Autoradiographic film

D. Related Products Available from Ambion

BrightStar™**Psoralen-Biotin**

Cat #1480

Nonisotopic labeling kit for fast, efficient biotin labeling of RNA, dsDNA, and oligonucleotides longer than 47-mers.

BrightStar™-Plus Membrane

Cat #10100–10104

Positively-charged nylon membrane recommended for use with Ambion's BrightStar™ nonisotopic labeling and detection products. These membranes are an excellent choice for Northern and other blot hybridizations.

II. BioDetect Protocol

A. When to Use BioDetect

Blot hybridizations (e.g. Northern, Southern, Dot Blots)

Biotinylated nucleic acids must be bound to positively-charged nylon membranes before beginning the detection procedure. This is typically achieved by hybridization of the probe with a target that is cross-linked to the blot.

Northern or Southern blots (preferably using BrightStar-Plus membrane from Ambion) on positively charged nylon membrane can be directly processed as described in section *II.C. Detection Procedure* on [page 4](#). Be sure to complete the post-hybridization washes before starting the BioDetect procedure. To control background, it is important to keep the membrane damp after the post-hybridization wash steps. Background may be higher using membranes that have been allowed to dry.

Gel separated probe (e.g. RPAs, SI Nuclease Assays)

The transfer of polyacrylamide gel-resolved probes to the BrightStar-Plus membrane, as in a ribonuclease protection assay, is the first step that differs from a typical ^{32}P -based assay. Electroblotting (instead of passive transfer) from the polyacrylamide gel onto the membrane is the best approach. After the transfer, the biotinylated nucleic acids are crosslinked to the membrane. Methods for both the membrane transfer and crosslinking are described in sections IV.A [on page 10](#), and IV.B [on page 11](#).

B. Buffer and Equipment Preparation

- The BioDetect protocol has been optimized for a 100 cm² membrane. Smaller or larger membranes can be used with appropriate (proportional) adjustments to solution volumes.
- The containers should be clean and RNase-free. We recommend treating the containers with Ambion's RNaseZap (Cat #9780), a strong RNase denaturing solution, followed by rinsing with high quality, RNase-free water.
- It is important that the Wash and Blocking Buffers are in solution for use. If any precipitate is visible, resolubilize it by warming to 37–65°C. Do not, however, heat the Streptavidin-Alkaline Phosphatase or the CDP-Star.

- Use good sterile technique when handling the Wash and Blocking Buffers. Bacterial contamination will cause high background.
- Dilute the needed amount of 5X Wash Buffer and 10X Assay Buffer to 1X with ddH₂O before use.
Prepare 5 ml of 1X Wash Buffer per cm² of membrane, or 500 ml of Wash Buffer for 100 cm² membrane.
Prepare 2 ml of 1X Assay Buffer per cm² of membrane, or 200 ml of Assay Buffer for 100 cm² membrane.
- Discard solutions after each incubation. The solutions are not reusable.

C. Detection Procedure



IMPORTANT!

All incubations described below should be done at room temperature with constant, gentle agitation in containers that allow free movement of the membrane (e.g. in flat trays).

Also, do not allow the membrane to dry! Once begun, the membrane must remain wet throughout the procedure.

I. Wash membrane 2 x 5 min in 1X Wash Buffer

Do two, 5 minute washes in approximately 1 ml Wash Buffer/cm² membrane/wash.

The Wash Buffer is provided at 5X, so it must be diluted prior to use. Generally, dilution in ddH₂O is sufficient (DEPC treatment and/or autoclaving of the water is not necessary).

The end of this washing period is good stopping point if there is not enough time to complete the detection procedure. At this point, the membrane can be stored wet, wrapped in plastic, for several days at 4°C.

2. Incubate membrane 2 x 5 min in Blocking Buffer

Do two, 5 minute incubations in approximately 0.5 ml Blocking Buffer/cm² membrane/incubation.

Do not dilute the Blocking Buffer as it is provided as a ready-to-use solution. As mentioned above, make sure to resuspend any precipitate prior to use.

3. Incubate membrane 30 min in Blocking Buffer


For this incubation, use approximately 1 ml Blocking Buffer/cm² membrane.

- 4. Incubate membrane
30 min in diluted
Strep-AP**

Prepare the diluted Strep-Alkaline Phosphatase (Strep-AP) by gently and thoroughly mixing together 10 ml Blocking Buffer and 1 μ l Strep-AP per typical size membrane (100 cm²).

Prepare only the volume needed for the experiment at hand, because Strep-AP is not stable for storage when it is diluted in Blocking Buffer.

It is important to mix the Strep-AP with the Blocking Buffer before adding it to the membrane, because if the Strep-AP is added directly to the container with the membrane and Blocking Buffer, heavy and uneven background may result.

 **NOTE:**
Vigorous agitation of enzymes can lead to partial denaturation.
- 5. Incubate membrane
10 min in Blocking
Buffer**

Use 0.5 ml Blocking Buffer/cm² membrane.

(optional) This incubation can be extended to 15 minutes to further decrease nonspecific background.
- 6. Wash membrane
3 x 5 min in 1X Wash
Buffer**

Do three, 5 minute washes in approximately 1 ml Wash Buffer/cm² membrane/wash.

(optional) These washes can be extended to 15 minutes each to further decrease nonspecific background.
- 7. Incubate membrane
2 x 2 min in 1X Assay
Buffer**

Do two, 2 minute incubations in approximately 0.5 ml 1X Assay Buffer/cm² membrane/incubation.

The Assay Buffer is provided at 10X, so it must be diluted in high quality water before use.
- 8. Incubate membrane
5 min in CDP-Star**

Use very little CDP-Star: only about 5 ml/100 cm². The CDP-Star is provided as a ready-to-use solution; do not dilute it.
- 9. Shake off excess
CDP-Star and expose
membrane to film**

Quickly blot the membrane on a piece of filter paper to remove excess CDP-Star, without letting the membrane dry. If any free-floating CDP-Star is left on the membrane, it may generate blotchy background. Wrap in a single layer of plastic, and expose to film at room temperature.
- 10. About the light
emission from
CDP-Star**

CDP-Star reaches peak light emission in 2–4 hours, then emission falls to a plateau which persists for several days. Try initial exposures of 1–2 hours followed by shorter exposures (five minutes to two hours) the following day if necessary to obtain the desired image.

III. Troubleshooting

A. Use of the Labeled Control DNA

The Psoralen-Biotin Labeled Control DNA is a linear 4 kb plasmid (1 ng/μl in 1X TE) containing a 1 kb insert of the mouse Glyceraldehyde-3-Phosphate Dehydrogenase (GAPDH) gene with biotin molecules covalently bound throughout. It is provided as a positive control for dot detections to ensure that the BioDetect Kit is functional. Typically, 100 fg to 1 pg of Labeled Control DNA can be detected using a dot detection assay.

Dot detection

To ensure the BioDetect Kit is performing well, detect dots of the Labeled Control DNA. To do this, make 10-fold serial dilutions of the Labeled Control DNA in molecular biology-grade water from 1 ng/μl to 100 fg/μl. Make the dots by simply pipetting 1 μl of each dilution onto dry BrightStar-Plus or another positively-charged nylon membrane. Allow the membrane to dry completely at room temperature.

Detect the dots by following the protocol described in section II.C. *Detection Procedure* starting on [page 4](#).

B. High Background

The most likely problems to occur in a nonisotopic detection protocol relate to background. Note that film exposure times using the BrightStar chemiluminescent system are significantly shorter (6 hours or less) than those when radiolabeled probes are used. An overnight exposure will produce high background. However if background comes up to unacceptable levels before the desired signal is detected, or the film is littered with spots and/or blotchiness, there are several possible sources and solutions for this problem that should be considered.

I. Precipitation of Wash or Blocking Buffers

The Wash and Blocking Buffers contain salts and detergents that may precipitate at room temperature. Accumulation of these precipitates on the membrane can result in increased background. Simply warming these solutions to between 37°C and 65°C to redissolve precipitates before use can prevent this problem.

Do *not* warm the Strep-AP or CDP-Star solutions.

- 2. Gel pieces and/or polyacrylamide residue on the membrane**

Directly after transfer, remove the membrane from the gel while submerged in transfer buffer. Then briefly rinse the membrane in the transfer buffer to wash off any loose gel fragments. Extending the length of washes in the detection procedure to fifteen minutes each (or longer) may also help.
- 3. Insufficient exposure to buffers and reagents**

If the membrane is not freely moving, it will not be fully exposed to the wash solutions, blocking buffers, and reagents. We have found that membranes processed in the conical tubes used in rotatory hybridization cabinets sometimes stick to the sides of these tubes. Although such tubes can be used successfully, we recommend using flat trays (e.g. the lids from pipet tip boxes). Washes following the incubation in diluted Strep-AP (steps [5–6](#) on [page 5](#)) may be extended to 15 minutes each to further decrease non-specific background when there is no other attributable source.
- 4. Failure to mix Strep-AP with Blocking Buffer before adding it to the membrane**

If the Strep-AP is added directly to the container with the membrane and Blocking Buffer, instead of mixing the solutions before pouring them into the container with the membrane, high background can result. This is probably due to localization of high concentrations of Strep-AP on the membrane surface. This problem can be avoided by thoroughly mixing the Strep-AP and Blocking Buffer before adding the solution to the membrane. We recommend mixing the Strep-AP and Blocking Buffer in a graduated cylinder with repeated, gentle inversion.
- 5. Excess CDP-Star not drained and blotted from membrane**

Since the CDP-Star reaction products are not bound by the alkaline-phosphatase, they can float freely on the membrane, creating blotchy background. After incubating the membrane with CDP-Star, hold the membrane up over the tray to allow any excess CDP-Star to drain off. Afterwards, briefly blot the membrane onto a piece of filter paper, making sure not to dry it.
- 6. Folds or multiple layers of plastic wrap**

Folds in the plastic wrap covering the membrane during exposure can trap CDP-Star causing “veins” and blotchiness. Also, multiple layers and plastic wrap edges can prevent the light from penetrating fully. Use a single plastic wrap layer and avoid creating wrinkles.
- 7. Insufficient liquid in filter papers during electroblotting**

The filter papers used to form the “sandwich” in the electroblotter must be dripping wet so that equal current can move through all parts of the transfer stack.

8. Bacterial/nuclease contamination

Poor signal detection coupled with an even, high background may indicate contamination of solutions and/or trays with nucleases or alkaline phosphatase. If you suspect contamination, thoroughly clean all vessels and trays with Ambion's RNaseZap (Cat # 9780) and rinse with RNase-free ddH₂O. Note, that if any of the solutions in the BioDetect Kit become contaminated with bacteria, they must be replaced.

C. Low Sensitivity

Low sensitivity usually results from problems preparing the probe or inefficient transfer of nucleic acids to the nylon membrane. The probe and the sensitivity of the detection procedure can be tested by spotting dilutions of the probe on a membrane strip, crosslinking, and following the detection protocol.

1. Low specific activity probe synthesized

At Ambion we get the highest specific activity biotinylated probes by post-synthesis labeling with Psoralen-Biotin. Ambion's BrightStar™ Psoralen-Biotin Labeling Kit (Cat #1480) can be used for this purpose.

The most likely factors that compromise specific activity of Psoralen-Biotin labeled probes are whether the correct wavelength was used for labeling and whether the UV source was too far from the sample. The purity of the sample and salt concentration can also be a factor.

Labeling reactions that incorporate biotinylated nucleotides enzymatically should contain a 2:3 ratio of biotin-CTP (or biotin-dCTP) to nonbiotinylated CTP or dCTP (a higher ratio will inhibit hybridization). Since biotinylated-NTPs are not incorporated as efficiently as unmodified nucleotides, further dilution will result in much less incorporation and a lower sensitivity probe. See Ambion's Technical Bulletin #173 for sources of modified NTPs and suggestions for setting up the reactions.

2. Inaccurate quantitation of probe

Contaminating salts, protein, and possibly acrylamide, can affect spectrophotometric readings. A_{260}/A_{280} ratios should fall between 1.8 and 2.1. If they do not, clean up biotinylated probes (e.g. by phenol extraction and/or precipitation and an EtOH wash step) to get a more accurate reading. Being aware of the maximum theoretical yield of a labeling reaction can also serve as a warning of inaccurate quantitation.

- 3. Degradation of probe** Probe degradation can result in lower sensitivity. Probe integrity can be assessed by running “probe alone” on a denaturing gel, transferring to membrane, crosslinking, and following the detection protocol.
- 4. Inefficient transfer to membrane** Inefficient transfer to membrane vs. “probe problems” can be assessed by staining the gel with EtBr after transfer to visualize any residual nucleic acid. In an RPA, if transfer is incomplete, a smear from the yeast RNA in the control lane should be visible; in a Northern, rRNAs should be visible. Alternatively, several transfers or similar dilutions of two probes transferred on the same membrane can be compared. Be sure to follow the manufacturer's instructions when electroblotting.
- 5. Inefficient or lack of crosslinking** Biotinylated probes transferred to a membrane from a nuclease protection assay gel must be crosslinked to the membrane before beginning the detection protocol; if left uncrosslinked, probes may wash off, reducing signal. However, this is not necessary for biotinylated probes used in blot hybridizations (e.g. Northern, Southern). These probes are bound by hybridization with the nucleic acid that is crosslinked to the membrane, and will not be removed by the BioDetect wash steps.
- 6. Exposure to film at reduced temperature** Exposure of the membrane to film below room temperature will inhibit or prevent the enzymatic reaction that produces light emission and will result in signal reduction or loss.

IV. Additional Procedures

A. Electroblothing from Polyacrylamide Gels

Electroblothing from polyacrylamide gels is best accomplished using a commercially available active transfer apparatus (tank electroblotter, semi-dry gel blotter, vacuum blotter, pressure blotter etc.). At Ambion we routinely use the OWL Scientific Panther™ Semi-Dry Electroblotter (Model # HEP-1), a relatively inexpensive and efficient electroblotter. The procedure is as follows:

1. Cut six pieces of filter paper and one BrightStar-Plus or other positively-charged nylon membrane to size of gel.
2. Thoroughly wet two pieces of the filter paper in 0.5X TBE and lay them over the bottom electrode. Separate the glass plates of the gel-plate sandwich and use a third piece of filter paper (wet) to lift the gel off the bottom plate.
3. Lay the wet membrane on the gel and wet the entire “sandwich” in 0.5X TBE, removing any bubbles that might have formed (a long plastic pipet can be rolled over the surface for this purpose).
4. Lay the “sandwich” carefully on the wet pieces of filter paper.
5. Thoroughly wet the remaining three pieces of filter paper and lay them carefully on the membrane surface. Again, it is important to avoid forming air bubbles.
6. Close the apparatus and apply current of 200 milliamperes for a period of 1 hour or 400 mA for 30 minutes.
7. After transfer is complete, remove the gel and membrane from the stack together and submerge them in transfer buffer. Peel the gel from the membrane and discard it. Rinse the membrane by briefly shaking it in the transfer buffer to dislodge any acrylamide fragments.
8. Immobilize the nucleic acid on the membrane by UV crosslinking (see below) and then proceed to the detection protocol (section *II.C. Detection Procedure* starting on [page 4.](#)).

B. Crosslinking

The membrane should be treated to immobilize the nucleic acid immediately after transfer and rinsing. This can be done by either one of the following methods:

UV crosslinking

If using a commercial crosslinker, the manufacturer's recommendations should be followed (i.e. UV exposure of approximately 1.2 millijoules). The membrane should be kept wet during this process. Ultraviolet crosslinking may also be accomplished by using a transilluminator or a handheld UV light source.

Baking

The membrane may be treated by baking at 80°C for 15 minutes. Since the membrane is not nitrocellulose, it is not necessary to use a vacuum oven. A conventional oven or convection oven is suitable.

V. Appendix

A. Specification Sheet

Kit Contents:

Amount	Component
1 L	5X Wash Buffer
3 L	1X Blocking Buffer
125 ml	10X Assay Buffer
20 μ l	Streptavidin-Alkaline Phosphatase
50 ml	CDP-Star®
20 μ l	Labeled Control DNA

Storage Conditions:

Store the 10X Assay Buffer and Streptavidin-Alkaline Phosphatase at 4°C.

Store the Labeled Control DNA at -20°C.

All other components should be stored at room temperature.

Quality Control:

The Labeled Control DNA is spotted onto Ambion's Bright-Star-Plus Membrane, and subjected to the procedure outlined in section *II.C. Detection Procedure* starting on [page 4](#) of this manual. Immediately after removing the CDP-Star, 10 pg of the Labeled Control DNA can be detected in a 30 min exposure to X-ray film. The membrane is exposed to x-ray film for 30 min again at 2 hr after removing the CDP-Star, and 1 pg of Labeled Control DNA can be detected.

Nuclease testing

Each component is tested in Ambion's rigorous nuclease assays.

Non-specific endonuclease/nickase activity

None detected after incubation with supercoiled plasmid DNA; analyzed on 1% agarose.

Exonuclease activity

None detected after incubation with ³²P-labeled *Sau3A* fragments of pUC19; analyzed by PAGE.

RNase activity

None detected after incubation with ³²P-labeled RNA; analyzed by PAGE.

Protease testing

None detected in protein-containing components after incubation with protease substrate, and analysis by spectrophotometry.

Safety:

This kit contains SDS and proprietary ingredients in the IX Blocking Buffer for which Material Safety Data Sheets are included on the following pages. The remaining kit components are dilute aqueous solutions containing buffers none of which are thought to present any health hazard.

B. IX Blocking Buffer Material Safety Data Sheet

Physical Data

Appearance and Odor	Liquid
Boiling Point	n/a
Specific Gravity	n/a
Vapor Density	n/a
Vapor Pressure (mm Hg)	n/a
Solubility in H ₂ O	Soluble

Fire and Explosion Hazard Data

Extinguishing Media	Water, CO ₂ , Foam, Chemical
Special Fire Fighting	Wear suitable protective gear for surrounding fire.
Fire/Explosion Hazards	May emit irritating and/or toxic fumes

Health Hazard Data

Effects of Overexposure	Irritant. May be harmful by inhalation or ingestion.
Emergency First Aid Procedures	Wash affected area with water. Irrigate eyes for at ≥15 min. See physician immediately. If ingested, drink several glasses of water, induce vomiting.

Reactivity Data

Stability	Stable
Incompatibility	Strong acids, bases or oxidants
Hazardous Decomp. Products	n/a
Hazardous Polymerization	n/a

Spill or Leak Procedures

If released or spilled	Ventilate area. Absorb spill with inert material. Clean area thoroughly.
Waste Disposal Method	Dispose according to federal, local and state regulations.

Special Protection Information

Respiratory Protection	Not expected to require personal respirator usage. (Use NIOSH approved respirator if necessary.)
Ventilation	Lab Hood for general ventilation
Protective Gloves	General lab safe gloves
Eye Protection	Use general eye protection-goggles
Handling and Storage	Wear appropriate protective clothing and gloves.

This bulletin is for your guidance and is based upon information and tests believed to be reliable. Ambion makes no guarantee of the accuracy or completeness of the data and shall not be liable for any damages thereto. The data are offered solely for your consideration, investigation, and verification. These suggestions should not be confused with either state, municipal, or insurance requirements, or with national safety codes and constitute no warranty. Any use of these data and information must be determined by the user to be in accordance with applicable federal, state, and local regulations.

C. SDS Material Safety Data Sheet

Physical Data

Appearance and Odor	Clear liquid, creates many bubbles when shaken.
Specific Gravity	0.35 g/cm ³
Solubility in H ₂ O	Soluble
CAS# 151-21-3	

Fire and Explosion Hazard Data

Flash Point	212°F
Extinguishing Media	Water, CO ₂ , Foam
Special Fire Fighting	Use self-contained breathing apparatus and protective clothing.
Unusual Fire/Explosion Hazards	May emit fumes of sulfur dioxide and carbon monoxide.

Health Hazard Data

Effects of Overexposure	May be harmful by inhalation, ingestion or skin absorption.
Threshold Limit Value	Unknown
Emergency First Aid Procedures	Flush with copious amounts of water. If ingested, notify physician and report that SDS has been ingested (detergent).

Reactivity Data

Stability	Stable
Incompatibility	Avoid heat and strong acids.
Hazardous Decomposition Products	Toxic fumes of carbon monoxide and sulfur dioxide
Hazardous Polymerization	Will not occur.

Spill or Leak Procedures

If released or spilled	Ventilate area. Place in containers with a lid. Wash spill area with copious amounts of water.
Waste Disposal Method	Review local and state regulations on appropriate disposal.

Special Protection Information

Respiratory Protection	Not expected to require personal respirator usage. (Use NIOSH approved respirator if necessary.)
Ventilation	Not expected to require any special ventilation.
Precautionary Labeling	Irritant
Handling and Storage Considerations	Use laboratory aprons and gloves. Store at room temperature.

This bulletin is for your guidance and is based upon information and tests believed to be reliable. Ambion makes no guarantee of the accuracy or completeness of the data and shall not be liable for any damages thereto. The data are offered solely for your consideration, investigation, and verification. These suggestions should not be confused with either state, municipal, or insurance requirements, or with national safety codes and constitute no warranty. Any use of these data and information must be determined by the user to be in accordance with applicable federal, state, and local regulations.